mikrobiologie labor technik

# Agar Chocolat

## PRINCIPIU

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
# Chocolate Agar

## PRINCIPLE

This medium is recommended for the growth of fastidious microorganisms such *Neisseria* and *Haemophilus*. Chocolate agar is adapted for the analysis of biological samples such as CRL and blood cultures.

## FORMULA

For preparation of chocolate agar can be used as a dehydrated medium: Columbia Agar Base or Blood Agar Base.

## METHOD

Melt the dehydrated medium, sterilize at 121°C, 15 minutes and cool to  $42 \pm 1^\circ\text{C}$ . Add 5 to 10% sheep or horse blood. Under slow and constant homogenization heat for 5 minutes at 70°C (to cook the blood). Cool for a second time and pour in sterile plates.

## PROCEDURE

Warm the plates to 37°C prior to inoculation. Incubate 24 to 48 hours under humid and possibly CO<sub>2</sub> enriched atmosphere.

## RESULTS

*Neisseria gonorrhoeae* colonies grow in 24 hours. They are small, greyish with regular contours and 0.5 to 1 mm Ø. If the incubation is prolonged these colonies grow and become irregular.

*Neisseria meningitidis* grow faster : they are slightly bigger with regular contours and with greyish and transparent coloration. Colonies can also be mucous.

*Neisseria* identification will be followed by a Gram coloration, oxydase test and a biochemical identification systems.

*Haemophilus* species grow easily on chocolate medium, specially when the atmosphere contain

CO<sub>2</sub>. Colonies are often mucous, translucent and greyish with a diameter of 0.5 to 1 mm.

## BIBLIOGRAPHY

1. Thayer J.D. and Martin J.E. 1966. Improved Medium for Cultivation of *N. gonorrhoeae* and *N. meningitidis*. Public Health Rep. 81:559-562.
2. Riou J.Y. et Guibourdenche M. 1992. Méthodes de laboratoires *Neisseria* et *Branhamella*. Inst. Past. – pp30-38.
3. COFRAC. Programme 116/00/BA24/00

## PACKAGING

### Columbia Agar Base dehydrated medium

Store between 1 and 30°C

140121A: Flask of 500 g

140121L: Flask of 500 g

### Blood Agar Base dehydrated medium

Store between 1 and 30°C

140131A: Flask of 500 g

140131L: Flask of 500 g

### Chocolate agar, ready to use plates

Store between 2 and 8°C

110176: Pack of 20 plates 90 mm Ø

120119: Pack of 32 tubes of 7 ml