mikrobiologie labor technik

Simmon's Citrat Agar

PRINCIPIU

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BIBLIOGRAFIE

1. Simmons J.S. 1926. A culture medium for differentiating organisms of typhoid-colon aerogenes groups and for isolating of certain fungi. *J. Infect. Dis.* 39:209-241.
2. Ewing W.H. and Edwards M.P.R. 1960. The principal divisions and groups of Enterobacteriaceae and their differentiation. *Int. Bull. Bact. Nomenclat. Taxon.* 10:1-12

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
Simmon's Citrate Agar

PRINCIPLE

Simmons Citrate Agar is used for differentiating Gram negative Bacilli through their ability to use citrate as a sole carbon source. The metabolism of citrate by the bacteria causes alkalinisation shown by the pH indicator turning the medium from green to blue, Bromthymol Blue.

FORMULA

Components	g/l
Magnesium sulfate	0.20
Ammonium dihydrogen phosphate	1.00
Dipotassium phosphate	1.00
Sodium citrate	2.00
Sodium chloride	5.00
Bromthymol Blue	0.08
Agar	15.00
Final pH : 6.9 ± 0.2 at 25°C	

METHOD

Suspend 24.2 grammes of dehydrated medium into one litre of purified water. Heat gently to boiling point under continuous stirring until completely dissolved. Dispatch 8 ml per tube. Autoclave for 15 minutes at 121°C. Cool in a slanted position with long slant and short butt.

PROCEDURE

With an inoculating needle or loop, pick the centre of a well isolated colony obtained from solid culture media. Streak only the surface of the slant with a light inoculum. Loosen the closure on the tube. Do not use a culture stemming from a broth since nutrients coming from the broth could

result in giving false positive. Incubate at 30 or 37°C for 1 to 7 days.

RESULTS

Observe daily the tubes. Citrate positive germs culture abundantly and turn the medium to blue. Citrate negative germs do not grow nor change the colour of the medium even after a long incubation period.

BIBLIOGRAPHY

1. Simmons J.S. 1926. A culture medium for differentiating organisms of typhoid-colon aerogenes groups and for isolating of certain fungi. J. Infect. Dis. 39:209-241.
2. Ewing W.H. and Edwards M.P.R. 1960. The principal divisions and groups of Enterobacteriaceae and their differentiation. Int. Bull. Bact. Nomenclat. Taxon. 10:1-12

PACKAGING

Dehydrated medium

(To be stored between 1 and 30°C)

140158A: Flask of 500 g

140158L: Flask of 500 g

Ready to use medium

110158: Pack with 20 plates of 90 mm Ø

120158: Pack with 5 tubes of 7 ml