mikrobiologie labor technik

T.S.I. Agar

PRINCIPIU

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
T.S.I. Agar

PRINCIPLE

T.S.I. (Triple Iron Sugar) is used for differentiating Gram-negative enteric bacilli based on the fermentation of dextrose, lactose and sucrose and on hydrogen sulfide production.

FORMULA

Components	g/l
Mix of peptone	20.00
Dextrose	1.00
Lactose	10.00
Sucrose	10.00
Sodium chloride	5.00
Sodium thiosulfate	0.20
Ferrous sulfate	0.20
Phenol red	0.025
Agar	13.00
Final pH: 7.3 + 0.2 at 25°C	

PREPARATION

Suspend 59,4 g in 1 liter purified water. Heat to boiling to dissolve completely. Dispense into tubes with closures. Autoclave at 118°C for 15 minutes. Cool in slanted position, as to have a slant and a butt of about 2 to 3 cm .

PROCEDURE

With an inoculating needle, prick the center of well-isolated colonies obtained from a solid media. Stab the center of the medium into the deep of the tube to 3 – 5 mm from the bottom. Withdraw the needle and streak the surface of the slant. Loosen closure on the tube before incubating 24 to 48 hours at 37°C. Read tubes for acid production of slant/butt, gas and hydrogen sulfide reactions.

RESULTS

- **An alkaline slant-acid butt (red/yellow)** indicates fermentation of dextrose only.

- **An acid slant-acid butt (yellow/yellow)** indicates fermentation of dextrose, lactose and/or sucrose.

- **An alkaline slant-alkaline butt (red/red)** indicates that neither dextrose nor lactose was fermented (non-fermentation).

- **Cracks, plits, or bubbles** in the medium indicate gas production.

- **A black precipitate** in the butt indicates hydrogen sulfide production.

REMARKUES

This medium, is similar to Kligler Iron Agar, it is added with sucrose that can be fermented faster than lactose by some gram-negative enteric bacilli. This allows to differentiate *Proteus* from *Salmonella*. It is common to sample a T.S.I. and a urea broth at the same time when wanting to identify *Proteus* or any other species that has an active urease. When T.S.I. has been prepared for more than a week before being used it is best to regenerate the medium by letting T.S.I. agar dissolve in hot water then cooling the tube as describe in the preparation paragraph.

LEGEND

AG : acid production (yellow) and gas

A : acid production (yellow)

NC : No colour change (red)

ALk : Alkaline reaction (red)

+ : Hydrogen sulfide production

- : No hydrogen sulfide production

	Butt	Slant	H2S
<i>Citrobacter</i>	AG	A	+
<i>Edwardsella</i>	A/AG	A	+
<i>Enterobacter aerogenes</i>	AG	A	-
<i>Enterobacter cloacae</i>	AG	A	-
<i>Escherichia coli</i>	AG	A	-
<i>Klebsiella</i>	AG	A	-
<i>Proteus mirabilis</i>	AG	A	+
<i>Proteus vulgaris</i>	AG	A	+
<i>Proteus morganii</i>	A/AG	NC/ALC	-
<i>Proteus rettgeri</i>	A	NC	-
<i>Shigella dysenteriae</i>	A	NC/ALC	-
<i>Shigella sonnei</i>	A	NC/ALC	-
<i>Salmonella typhi</i>	A	NC/ALC	+
<i>Salmonella paratyphi A</i>	AG	NC/ALC	-
<i>Salmonella choleraesuis</i>	AG	NC/ALC	-
<i>Salmonella enteritidis</i>	AG	NC/ALC	+
<i>Salmonella typhimurium</i>	AG	NC/ALC	+

BIBLIOGRAPHY

1. American Public Health Association. 1963. Diagnostic Procedures and Reagents. 4th Ed. pp. 150 and 294-295. A.P.H.A. Inc., New York.
2. American Public Health Association. 1966. Recommended Methods for the Microbiological Examination of Foods. 2nd Ed. pp. 158 and 185, A.P.H.A. Inc., New York.
3. Hajna A.A. 1945. Triple-Sugar Iron Medium for the Identification of the Intestinal Group of Bacteria. J. Bact., 49:516-517.
4. US Pharmacopoeia. XVI medium.
5. European Pharmacopoeia. M medium.

PACKAGING

Dehydrated medium

(Store between 1 -30°C)

1401120A: Flask of 500 g

1401120L: Flask of 500 g

Prepared medium in flask

(Store between 2 25°C)

120117: Pack of 32 slanted tubes