mikrobiologie labor-technik

Müller Kauffman Bulion

PRINCIPIU

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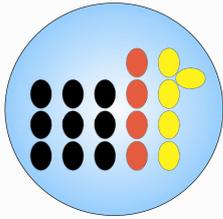
Müller Kauffman Broth

PRINCIPLE

Müller-Kauffmann broth's formula corresponds to the one described by Müller, and then modified by Kauffmann. This fluid medium is used as a selective enrichment broth when screening for *Salmonella* in foodstuffs, water and biological specimens.

The medium is composed of sodium thiosulfate that is oxidised into tetrathionate by a solution of iodine/iodide added to the base at the last minute. This contributes to inhibit the growth of coliforms and most of all bacteria belonging to the intestinal flora. *Proteus* and *Salmonella* can reduce tetrathionate and therefore find their intake of sulfur for their growth. Calcium neutralises the formed sulfuric acid in order to prevent any drop of pH levels that might compromise the growth of bacteria. Brilliant green inhibits the growth of gram positive. Oxgall contributes to the growth of *Salmonella* by slowing down the growth of interfering flora.

FORMULA

Components	g/l
Tryptone	7.00
Soy peptone	2.30
Sodium chloride	2.30
Calcium carbonate	25.00
Sodium thiosulfate 5H ₂ O	40.70
Oxgall	4.75

Iodine/iodide Solution

Iodine	20.00 g
Potassium Iodide	25.00 g
Purified sterile water	100.00 ml

Dissolve the potassium Iodide in 5 ml of purified water, add iodine and heat slightly

to help to dissolve. Complete to 100 ml with purified water.

Brilliant Green 0.1% solution.

Brilliant green	0.10 g
Purified water	100.00 ml

Suspend brilliant green powder in to the purified water, homogenise. Heat 30 minutes at 100°C under constant stirring. Keep the prepared solution in a brown flask.

PREPARATION

Suspend 82 g of powder in 1 litre of purified water. Boil for one minute, under continuous stirring until completely dissolved. Cool to 45-50°C. Add 19 ml of Iodine/Iodide solution and 9.5 ml of 0.1 % Brilliant green solution. Homogenise well and dispatch 10 ml per tube.

PROCEDURE

Inoculate 10 ml (or 20 ml) tubes with 1 ml (or 2 ml) of sample primarily enriched in non selective broth. Homogenise well then incubate at 41-43°C for 18-24 hours or up to 48 hours depending on the procedure carried out.

RESULTS

After incubation periods, subculture onto an adequate selective agar (ASAP, Hektoën, X.L.D . or modified XLT4).

LIMITS AND PRECAUTIONS

In order to keep active the selective properties of the complete medium Brilliant green solution must be added to the Müller-Kauffmann base according to the preparation specifications. Müller-Kauffmann broth contains carbonate calcium that is used as a buffering agent which tends to sediment at the bottom of

tubes and flasks. Tubes and flasks have to be homogenised before inoculation. Complete medium has a shelf life of one week. It is wise to use a second selective broth in parallel to Müller-Kauffmann such as selenite cystine which would be incubated at 37°C or Rappaport Vassiliadis incubated at 42°C.

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